

This kit is intended for Research Use Only.

Not for use in diagnostic procedures.

NAME AND INTENDED USE

Anti-Parietal Cell is an indirect solid phase enzyme immunoassay (ELISA) for measurement of IgG class autoantibodies against a- and b-subunits of the Parietal Cell H⁺/K⁺-ATPase in human serum or plasma.

PRINCIPLE OF THE TEST

Highly purified pig Parietal Cell H⁺/K⁺-ATPase is bound to microwells. Antibodies against this antigen, if present in diluted serum or plasma, bind to the respective antigen. Washing of the microwells removes unspecific serum and plasma components. Horseradish peroxidase (HRP) conjugated anti-human IgG immunologically detects the bound specimen sample antibodies forming a conjugate/antibody/antigen complex. Washing of the microwells removes unbound conjugate. An enzyme substrate in the presence of bound conjugate hydrolyzes to form a blue color. The addition of an acid stops the reaction forming a yellow end-product. The intensity of this yellow color is measured photometrically at 450 nm. The amount of colour is directly proportional to the concentration of IgG antibodies present in the original sample.

WARNINGS AND PRECAUTIONS

1. All reagents of this kit are strictly intended for Research Use Only.
2. Do not interchange kit components from different lots.
3. Components containing human serum were tested and found negative for HBsAg, HCV, HIV1 and HIV2 by FDA approved methods. No test can guarantee the absence of HBsAg, HCV, HIV1 or HIV2, and so all human serum based reagents in this kit must be handled as though capable of transmitting infection.
4. Avoid contact with the TMB (3,3',5,5'-Tetramethylbenzidine). If TMB comes into contact with skin, wash thoroughly with water and soap.
5. Avoid contact with the Stop Solution which is hydrochloric acid (1 M). If it comes into contact with skin, wash thoroughly with water and seek medical attention.
6. Some kit components (i.e. Controls, Sample buffer and Buffered Wash Solution) contain Sodium Azide as preservative. Sodium Azide (NaN₃) is highly toxic and reactive in pure form. At the product concentrations, though not hazardous. Despite the classification as non-hazardous, we strongly recommend using prudent laboratory practices (see 8., 9., 10.).
7. Some kit components contain Proclin 300 as preservative. When disposing reagents containing Proclin 300, flush drains with copious amounts of water to dilute the components below active levels.
8. Wear disposable gloves while handling specimens or kit reagents and wash hands thoroughly afterwards.
9. Do not pipette by mouth.

DRG® Anti-Parietal Cell (EIA-3607)



Revised 11 Jan. 2012 rm (Vers. 1.1)

RUO	in the USA
-----	------------

10. Do not eat, drink, smoke or apply makeup in areas where specimens or kit reagents are handled.
11. Avoid contact between the buffered Peroxide Solution and easily oxidized materials; extreme temperature may initiate spontaneous combustion.

Observe the guidelines for performing quality control in medical laboratories by assaying controls and/or pooled sera. During handling of all kit reagents, controls and serum samples observe the existing legal regulations.

CONTENTS OF THE KIT

Package size	96 determ.
Qty.1	Divisible microplate consisting of 12 modules of 8 wells each, coated with highly purified pig H ⁺ /K ⁺ -ATPase, a- and b-subunits. Ready to use.
6 vials, 1.5 ml each	Standards: Anti-Parietal Cell Calibrators (A-F) in a serum/buffer matrix (PBS, BSA, NaN ₃ <0,1% (w/w)) containing: 0; 6.3; 12.5; 25; 50; 100 U/ml. Ready to use.
2 vials, 1,5 ml each	Anti-Parietal Cell Controls in a serum/buffer matrix (PBS, BSA, NaN ₃ <0,1% (w/w)) positive (1) and negative (2), for the respective concentrations see the enclosed package insert. Ready to use.
1 vial, 20 ml	Sample buffer (Tris, NaN ₃ <0,1% (w/w)), yellow, concentrate (5x).
1 vial, 15 ml	Enzyme conjugate solution (PBS, PROCLIN 300 <0,5% (v/v)), (light red) containing polyclonal rabbit anti-human IgG; labelled with horseradish peroxidase. Ready to use.
1 vial, 15 ml	TMB substrate solution. Ready to use.
1 vial, 15 ml	Stop solution (1 M hydrochloric acid). Ready to use.
1 vial, 20 ml	Wash solution (PBS, NaN ₃ <0,1% (w/w)), concentrate (50x).

STORAGE AND STABILITY

1. Store the kit at 2-8 °C
2. Keep microplate wells sealed in a dry bag with desiccants
3. The reagents are stable until expiration of the kit
4. Do not expose test reagents to heat, sun or strong light during storage and usage
5. Diluted sample buffer and wash buffer are stable for at least 30 days when stored at 2-8 °C

MATERIALS REQUIRED**Equipment**

- Microplate reader capable of endpoint measurements at 450 nm
- Multi-Channel Dispenser or repeatable pipet for 100 µl
- Vortex mixer
- Pipets for 10 µl, 100 µl and 1000 µl
- Laboratory timing device
- data reduction software

Preparation of reagents

- distilled or deionized water
- graduated cylinder for 100 and 1000 ml
- plastic container for storage of the wash solution

SPECIMEN COLLECTION, STORAGE AND HANDLING

1. Collect whole blood specimens using acceptable medical techniques to avoid hemolysis
2. Allow blood to clot and separate the serum by centrifugation
3. Test serum should be clear and non-hemolyzed. Contamination by hemolysis or lipemia is best avoided, but does not interfere with this assay.
4. Specimens may be refrigerated at 2-8 °C for up to five days or stored at –20 °C up to six months.
5. Avoid repetitive freezing and thawing of serum samples. This may result in variable loss of autoantibody activity
6. Testing of heat-inactivated sera is not recommended

PROCEDURAL NOTES

1. Do not use kit components beyond their expiration dates
2. Do not interchange kit components from different lots
3. All materials must be at room temperature (20-28 °C)
4. Have all reagents and samples ready before start of the assay. Once started, the test must be performed without interruption to get the most reliable and consistent results.
5. Perform the assay steps only in the order indicated
6. Always use fresh sample dilutions
7. Pipette all reagents and samples into the bottom of the wells
8. To avoid carryover contamination change the tip between samples and different kit controls
9. It is important to wash microwells thoroughly and remove the last droplets of wash buffer to achieve best results.
10. All incubation steps must be accurately timed
11. Control sera or pools should routinely be assayed as unknowns to check performance of the reagents and the assay.
12. Do not re-use microplate wells

For all controls, the respective concentrations are provided on the labels of each vial. Using these concentrations a calibration curve may be calculated to read off the specimen sample results semi-quantitatively.



Revised 11 Jan. 2012 rm (Vers. 1.1)

RUO in the USA

PREPARATION OF REAGENTS

Preparation of sample buffer

Dilute the contents of each vial of the sample buffer concentrate (5x) with distilled or deionized water to a final volume of 100 ml prior to use. Store refrigerated: stable at 2-8 °C for at least 30 days after preparation or until the expiration date printed on the label.

Preparation of wash solution

Dilute the contents of each vial of the buffered wash solution concentrate (50x) with distilled or deionized water to a final volume of 1000 ml prior to use. Store refrigerated: stable at 2-8 °C for at least 30 days after preparation or until the expiration date printed on the label.

Sample preparation

Dilute all specimen samples **1:100** with sample buffer before assay. Therefore combine 10 µl of sample with 990 µl of sample buffer in a polystyrene tube. Mix well. Controls are ready to use and need not be diluted.

TEST PROCEDURE

1. Prepare a sufficient number of microplate modules to accommodate controls and prediluted specimen samples.
2. Pipet 100 µl of calibrators, controls and prediluted specimen samples in duplicate into the wells.

	1	2	3	4	5	6
A	SA	SE	P1	P5		
B	SA	SE	P1	P5		
C	SB	SF	P2	P..		
D	SB	SF	P2	P..		
E	SC	C1	P3			
F	SC	C1	P3			
G	SD	C2	P4			
H	SD	C2	P4			

SA-SF: standards A to F
 P1, P2...C: specimen sample 1, 2 ...
 C1: positive control
 C2: negative control

3. Incubate for 30 minutes at room temperature (20-28 °C)
4. Discard the contents of the microwells and wash 3 times with **300 µl** of wash solution.
5. Dispense **100 µl** of enzyme conjugate into each well
6. Incubate for 15 minutes at room temperature
7. Discard the contents of the microwells and wash 3 times with **300 µl** of wash solution
8. Dispense **100 µl** of TMB substrate solution into each well
9. Incubate for 15 minutes at room temperature
10. Add **100 µl** of stop solution to each well of the modules and incubate for 5 minutes at room temperature
11. Read the optical density at 450 nm and calculate the results. Bi-chromatic measurement with a reference at 600-690 nm is recommended.

DRG[®] Anti-Parietal Cell (EIA-3607)

Revised 11 Jan. 2012 rm (Vers. 1.1)

RUO	in the USA
-----	------------

The developed colour is stable for at least 30 minutes. Read optical densities during this time.

Automation

The Anti-Parietal Cell ELISA is suitable for use on open automated ELISA processors. The test procedure detailed above is appropriate for use with or without automation.

Calculation of results

For Anti-Parietal Cell a 4-Parameter-Fit with lin-log coordinates for optical density and concentration is the data reduction method of choice.

Recommended Lin-Log Plot

First calculate the averaged optical densities for each calibrator well. Use lin-log graph paper and plot the averaged optical density of each calibrator versus the concentration. Draw the best fitting curve approximating the path of all calibrator points. The calibrator points may also be connected with straight line segments. The concentration of unknowns may then be estimated from the calibration curve by interpolation.

Calculation example

The figures below show typical results for Anti-Cardiolipin IgG/IgM ELISA. These data are intended for illustration only and should not be used to calculate results from another run.

Calibrators									
No	Position	OD 1	OD 2	Mean	Conc. 1	Conc. 2	Mean	decl. Conc.	CV %
STA	A 1/B 1	0.016	0.016	0.016	0.00	0.00	0.00	0.0	0.0
STB	C 1/D 1	0.332	0.335	0.334	6.45	6.40	6.43	6.3	0.6
STC	E 1/F 1	0.548	0.558	0.553	12.0	12.2	12.1	12.5	1.3
STD	G 1/H 1	0.934	0.956	0.945	25.6	24.8	25.2	25.0	1.6
STE	A 2/B 2	1.410	1.386	1.398	51.0	50.0	50.5	50.0	1.2
STF	C 2/D 2	1.823	1.840	1.832	98.3	101.1	99.7	100.0	1.8

REFERENCES

1. Irvine, W. J., Davies, S. H., Delamore, I. W., Williams, A. W.. Immunological relationship between pernicious anemia and thyroid disease.
Br. Med. J. 1962, pp. 454-456.
2. Taylor, K. B., Roitt, I. M., Doniach, D., Couchman, K. G., Shapland, C.. Autoimmune phenomena in pernicious anemia: gastric antibodies.
Br. Med. J. 1962, pp. 1347-1352.
3. Rabon, E. C. and Reuben M. A.. The mechanism and structure of the gastric H/K-ATPase.
Annu. Rev. Physiol. 1990, No. 52, pp. 321-344.
4. Pedersen, P. L. and Carfoli, E.. Ion motive ATPases. I. Ubiquity properties and significance to cell function.
Trends Biochem Sci 1987, No. 12, pp. 146-149.
5. Van Driel, I. R., Callaghan, J. M.. Proton and potassium transport by H/K ATPases.
Clin. Exp. Pharmacol. Physiol. 1995.
6. Irvine, W. J., Cullen, D. R., Mawhinney, H.. Natural history of autoimmune achlorhydric atrophic gastritis. A 1-15 year follow up study.
Lancet 1974, No. 2, pp. 482-485.
7. Whittingham, S., Mackay, I. R.. Pernicious anemia and gastric atrophy. In: Rose, N. R., Mackay, I. R. eds. The Autoimmune Diseases. New York, Academic Press 1985, pp. 243-266.
8. Uibo, R., Krohn, K., Villako, K., Tammur, T., Tamm, A.. The relationship of parietal cell, gastrin cell and thyroid autoantibodies to the state of the gastric mucosa in a population sample. Scand.
J. Gastroenterol. 1984, No. 19, pp. 1075-1080.
9. Varis, K., Ihamako, T., Harkonen, M., Samloff, I. M., Siurala, M.. Gastric morphology, function and immunology in first degree relatives of probands with pernicious anemia and controls.
Scand. J. Gastroenterol. 1979, No. 14, pp. 129-139.